

# pH-Fix 0-14

en

## Test strips for the determination of the pH-Value in the range of pH 0-14

### Contents:

- 100 Test strips
- 1 Colour scale pH 0-14

### Instructions for use:

1. Dip test strip for 1 sec. with all test fields in the sample solution.
2. Shake off excess sample solution.
3. Compare with colour scale and read off the corresponding pH-value.

### Note:

If sample solution is weakly buffered leave test strip in solution until colour no longer changes.

### Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container below + 30 °C in a dry place.

### Disposal:

Please dispose used test strips in domestic waste.

